STERYFLON PLUS

PTFE membrane cartridge



- High permeability versus gas
- Thermic cycles resistant
- Repeatedly steamable in situ or in autoclave
- Thermowelded construction
- FDA-listed materials per CFR21
- **Bio-Safety per USP-Plastics**
- Validation Guide available on request

STERYFLON Plus is a filter element designed, manufactured and submitted to severe controls for applications in processes where the effluent is required to be free from particellar and biological contaminants.

(EN) Data sheet

The expanded PTFE membrane, intrinsically hydrophobic, allows superior performances both in gas filtration and non-acqueos liquid filtration. Bio-tech pharmaceutical, electronic, food & beverage industries can roly on a product with high standard of quality suitable to solve contamination issues on critical applications.

STERYFLON is manufactured within a controlled environment and each cartridge is integrity tested and is validated for bacterial retention correlated with microbiological challenge test.

The retention is checked on regular sampling.

MATERIALS OF CONSTRUCTION

Filter media	PTFE membrane	
Upstream supports	orts polypropylene	
Downstream supports	polypropylene	
Internal Core	polypropylene	
External Cage	polypropylene	
End caps / Adapters	polypropylene	

FOOD-SAFETY

STERYFLON Plus filter elements meet regulation (EC) 1935/2004 for indirect food contact. **BIO-SAFETY**

Filter media and components pass USP Biological Reactivity and Chemical-Physical tests for CLASS VI plastics. The filter meets USP "Water for injection" requirements for particle release and the effluent is Non-Reactivity

particle release and the effluent is Non-Pyrogenic per USP Bacterial Endotoxins (< 0,25 EU/ml). QUALITY STANDARDS

Produced under a certified Quality System to guarantee traceability of manufacturing records and integrity testing results.

OPERATING CONDITIONS

- max. continuous temperature of fermentation inlet and exhaust air	70 °C
- max. continuous temperature for vent filter in recirculation loop	83 °C
- max. cumulative time of steam sterilization	150 hours at 140 °C with cycles of 30 minutes
- sanitization with chemicals	can be sanitized by standard chemical agents
- max. differential pressure	5,0 bar at 25 °C
- recommended change out differential pressure	2,0 bar at 25 °C

CODE	ABSOLUTE FILTRATION RATING		3) BACTERIAL RETENTION >10 ⁷ CFU/cm ²	ACCEPTABLE LIMIT FOR INTEGRITY TEST
	IN LIQUID	IN DRY GAS		FOR 10" CARTRIDGE
SH	0,1 µm	< 0,01 µm	Acheleoplasma laidlawii in liquid	1) ≤ 25 ml/min @ 1,0 bar
SL	0,2 µm	< 0,01 µm	B. Diminuta in liquid / Batteriofago T1 in aerosol	1) ≤ 12 ml/min @ 0,8 bar
SLA	0,2 µm	< 0,01 µm	Brevundimonas diminuta in liquid	2) ≤ 16 Nml/10min @ 2,5 bar
NI J			-	·

Note:

sale

1- The integrity is verified by DIFFUSION TEST using an aqueous solution of isopropyl alcohol (IPA 60/40 V/V) as wetting liquid.
2. The integrity is used to be a solution of the integrity of the integrity of the integrity.

2- The integrity is verified by WATER FLOW INTRUSION TEST.

comm

3- Cartridges are validated for retention of the microorganism reported in the table according to ASTM F838.











We are here to help. Contact us!






STERYFLON PLUS

PTFE membrane cartridge


EN Data sheet



AIR FLOW RATE FOR 10" CARTRIDGE

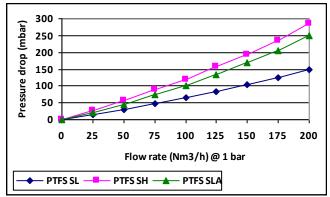

PTFS - <u>207</u> <u>1</u>

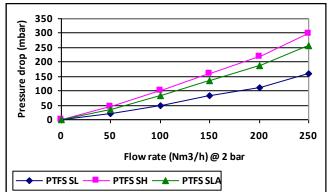

STERYFLON Plus ORDERING INFORMATION

END FITTING	CODE
DOE: double open end with flat gaskets	200
SOE: open end with (2) O-Ring 2.222. Blind end with flat top.	203
SOE: open end with (2) O-Ring 2.226 and 2 bayonet locks. Blind end with fin.	207
SOE: open end with (2) O-Ring 2.222. Blind end with fin.	208
SOE: open end with (2) O-Ring 2.222 and 3 bayonet locks. Blind end with fin.	212

ABSOLUTE FILTRATION RATING micron	CODE
0,1	SH *
0,2	SL *
0,2	SLA **

<u>SL</u>

-

* Integrity by IPA diffusion test ** Integrity by water intrusion test

CODE NOMINAL
LENGTH
05 5″
1 10"
2 20"
3 30″
4 40″

CODE	PRODUCT GRADE	
No code	Biological Grade	
рн	Non-pyrogenic Grade. Quality Certification in the box	
рнн	Non-pyrogenic Grade. Quality Certification, with serial number, in the box	

<u>PH</u>

CODE	GASKETS		END FITTING	
No code	Standard	EPDM	200	
т	On request	Teflon	200	
No code	Standard	Silicone		
v	On request	Viton	All the others	
E	On request	EPDM		
F	On request	FEP	207	







